


Experience convenience and quality while reducing your lab's waste by 30%

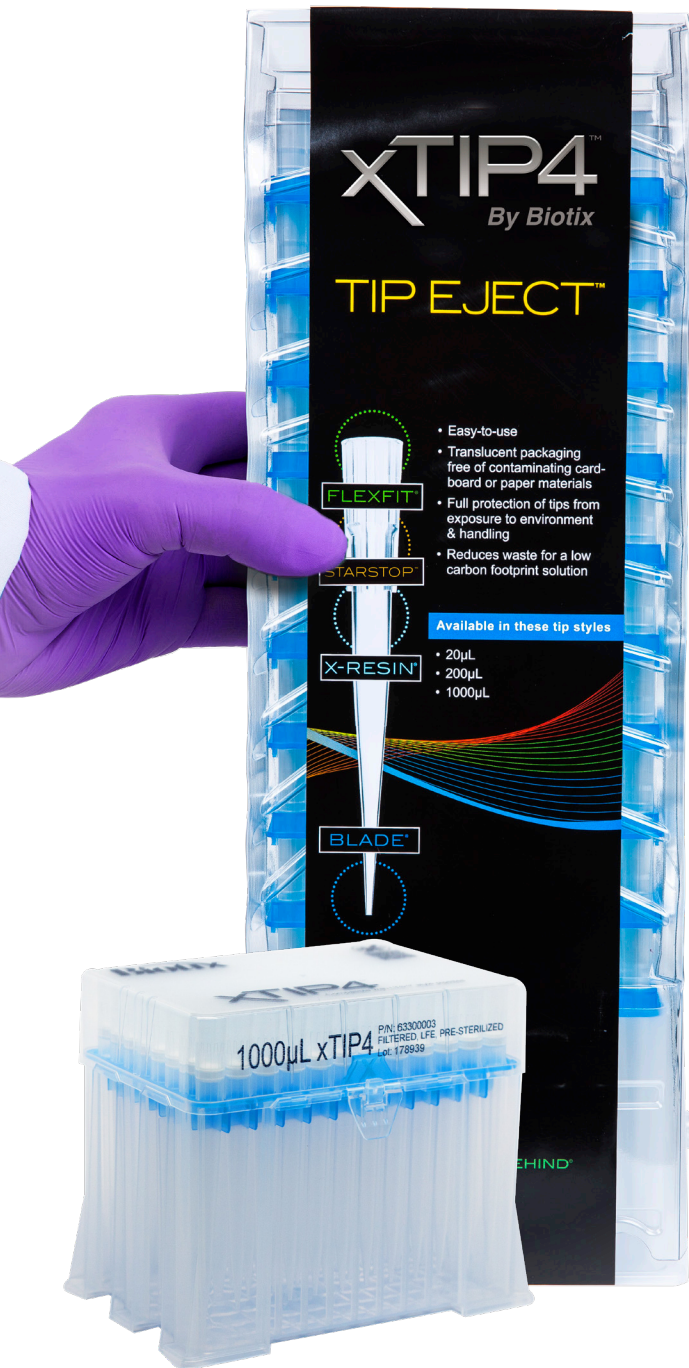


This self-aligning reload solution allows you to easily reload your racks without coming into contact with the tips. The rigid plastic covering also shields unused tips from contamination for later use.

- Available only for LTS and xPIPETTE compatible tips.
- Vertically stacked snapcards take up less space on your lab bench.
- Reduces plastic waste by 30% when compared to racked tips.

Pipette tip racks are fully autoclaveable at 121° C after transfer from reload.



Learn more about our sustainability initiatives at [BIOTIX.com](http://BIOTIX.com)

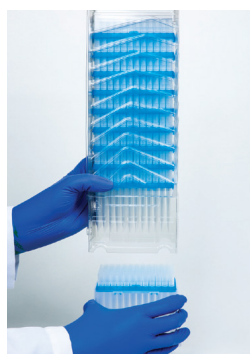


**Step 1: Remove the base**

Remove the lower protective rack base.



**Step 2: Align empty rack**

Align and place Tip Eject over your empty Biotix rack.

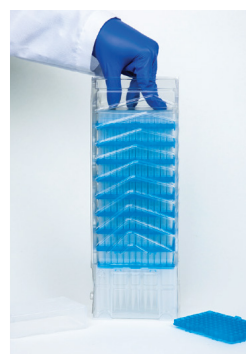

**Step 3: Push down**

Push down until a “click” is heard, indicating tips are locked into rack.



**Step 4: Replace stack over empty rack base**

Repeat to refill empty racks as needed

Volume $\mu\text{L}$	Pre-Sterilized	xTIP Tip Eject Reload	xTIP Tip Eject Packaging
10	Yes		
	No		
10 XL	Yes		
	No		
20	Yes	12111386	10 refills/pack
	No		
200	Yes		
	No		
250	Yes	12111387	10 refills/pack
	No		
300	Yes		
	No		
1000	Yes	12111389	8 refills/pack
	No		
1200	Yes	12111390	8 refills/pack
	No		

Distributed by Fisher Scientific. Contact us today:

**In the United States**

Order online: [fishersci.com](http://fishersci.com)

Call customer service: 1-800-766-7000

**In Canada**

Order online: [fishersci.ca](http://fishersci.ca)

Call customer service: 1-800-234-7437

